

QUICK REFERENCE GUIDE

Record the MAX & MIN temperatures twice daily - early morning & late afternoon. This will show what has happened during the day & overnight:

Recording the MINIMUM & MAXIMUM temperatures

- Press the 'I' button once to display 'Thi' (Temp High)
- Then press and hold "i" again to display MAX temperature
- Record the displayed MAX temperature
- Release button to display 'Tlo' (Temp Low)
- Then press and hold "i" again to display MIN temperature
- Record the displayed MIN temperature, then release "i".
- After this 'T1' will be displayed – ignore this. The current temperature will be displayed after a few seconds

Resetting the MINIMUM & MAXIMUM temperatures

- *Repeat above then :*
- When pressing & holding "i" to display the MAX temperature press 'X' at the same time
- This has now cleared the MAX temperature memory and reset it to the current temperature
- When pressing & holding "i" to display the MIN temperature press 'X' at the same time
- This has now cleared the MIN temperature memory and reset it to the current temperature
- After this 'T1' will be displayed – ignore this. The current temperature will be displayed after a few seconds

If you do not press either buttons for a few seconds the controller will revert back to displaying the internal temperature as normal.

Alarm Display

If the temperature falls below +2c or rises above +8c for more than 15 minutes, the alarm will be activated. 'Hi' or 'Lo' will flash in the display and the buzzer will sound. Press the 'X' button to silence the buzzer.

Automatic Defrost

During the automatic defrost period 'def' will displayed. Normal storage temperatures are maintained during this defrosting period.